
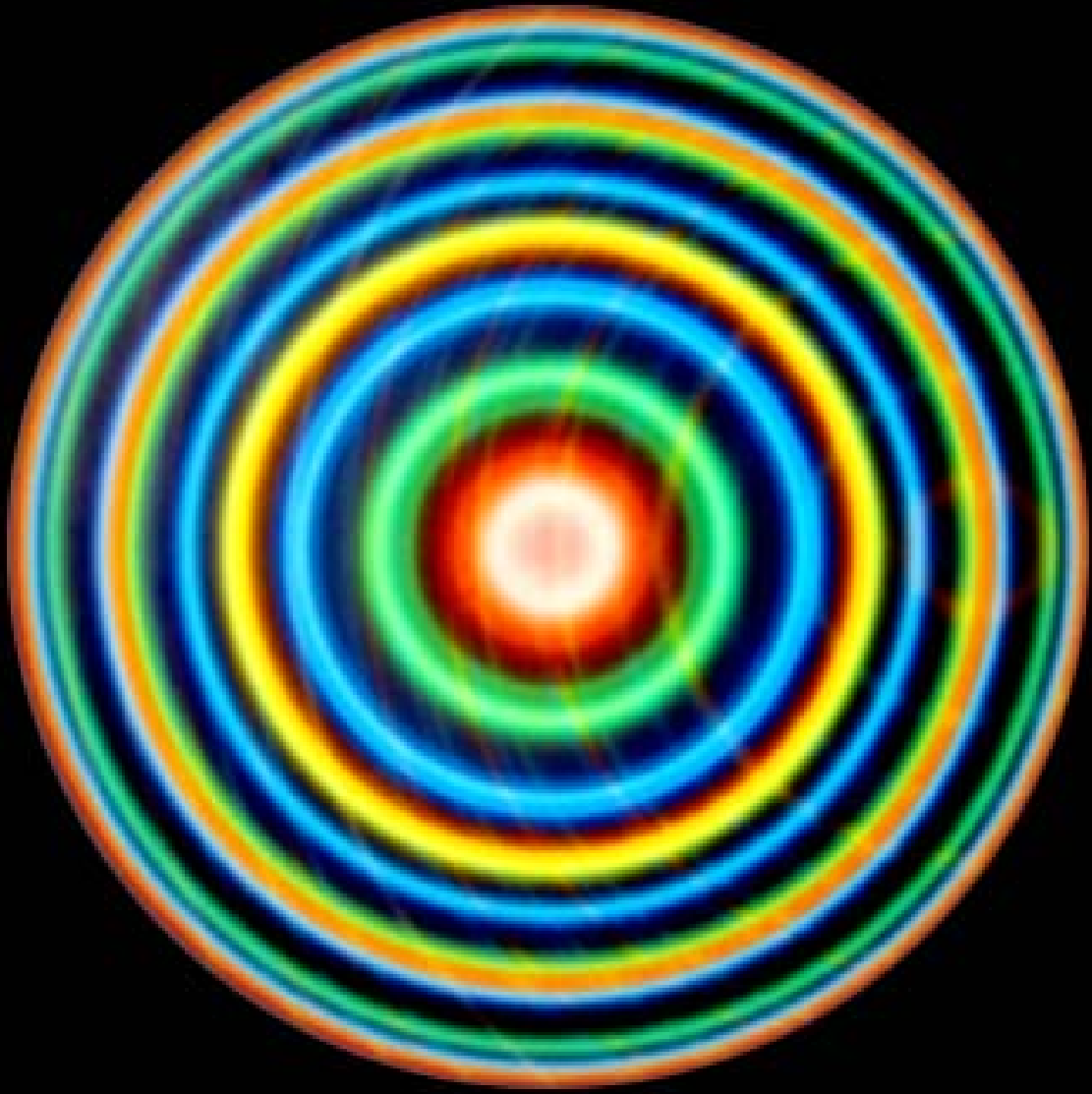
Prismen

Bernhard Halle Nachfl.

90°-Prismen

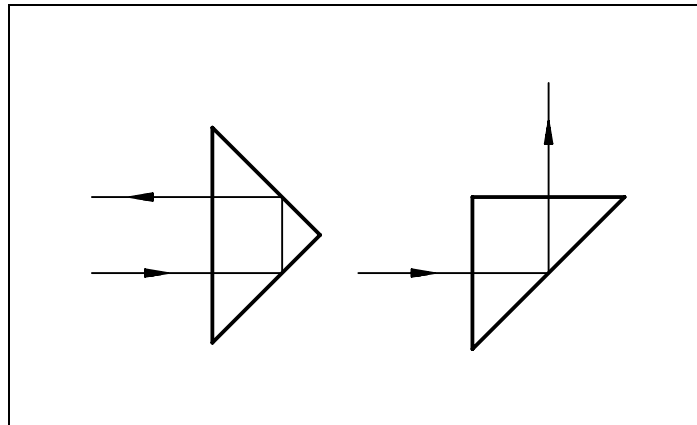Right Angle Prisms

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The prisms are made of crown glass BK 7 or fused silica Suprasil 1.

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Quality B:

Angular error up to approx. 2'. Flatness of surfaces up to $\lambda/5$. All edges beveled.

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
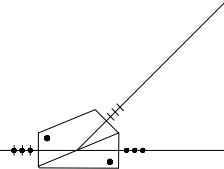
Antireflection and other coatings on request
Other materials and sizes on request

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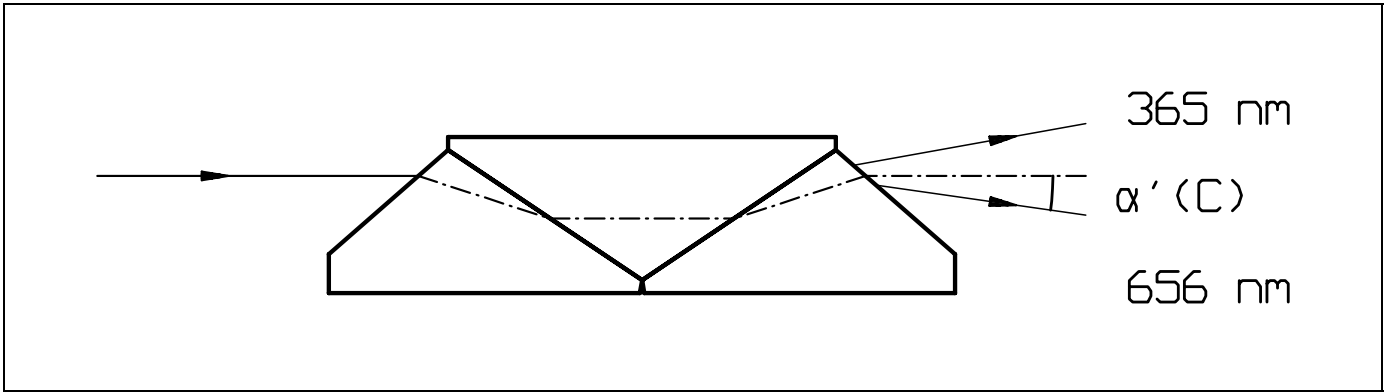This direct vision dispersing prism consists of three optically contacted parts. The outer prism parts are fused silica Suprasil, whereas the material in the inner prism is a flint glass with high UV transmission. The internal transmittance is 80% for a 20 mm glass path at the wavelength 365 nm. Thus in lower light power situations the prism can be used up to the near UV region (above about 350 nm).

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The direct vision is at the wavelength $\lambda = 436$ nm, the angular dispersion $\alpha_F - \alpha_C$ is $4,9^\circ$.

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Standard version:

The usable beam cross-section is about 6×10 mm² for a prism with a cross-section of 18×10 mm² and a length of 72 mm. DAM 610

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The Amici prism can also be supplied mounted. additional price DAM 0

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Available on special order:

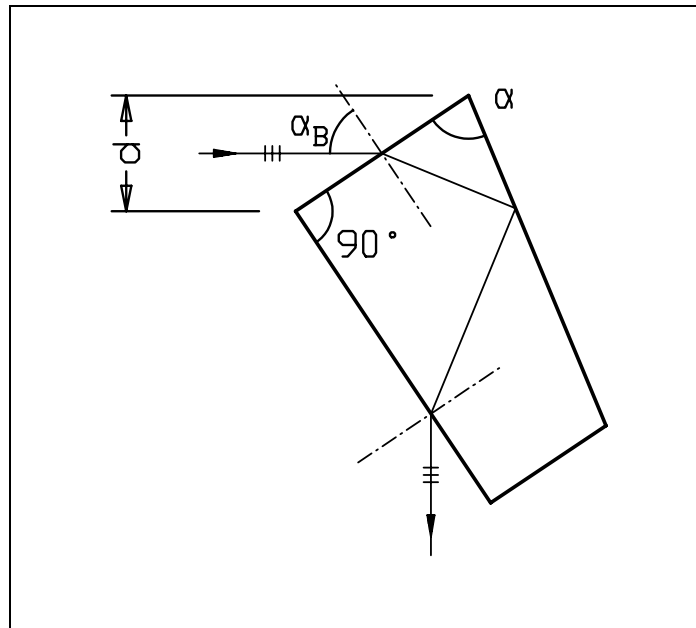
Two-part Amici prisms with direct vision wavelength and dispersion according to your specification on request

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more-part Amici prisms with direct vision wavelength and dispersion according to your specification on request

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The Brewster dispersion prisms are used for dispersion and 90° deviation of linearly polarized light (p-polarization), with no reflection losses. It consists of one solid glass block with three polished surfaces.

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The angle α is calculated so that the wave polarized parallel to the plane of incidence (p-polarization) completely enters the prism and then exits it again without any reflection losses, when the total deviation is 90° . This condition can only be achieved accurately for one wavelength.

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
The three polished faces are plane to less than $\lambda/10$. The angle errors are less than $2'$.

δ	R_p (280 nm)		R_p (200 nm)	
	$\alpha_B + \delta$	$\alpha_B - \delta$	$\alpha_B + \delta$	$\alpha_B - \delta$
0.00 °	0 %	0 %	0.01114 %	0.01114 %
0.05 °	0.00004 %	0.00002 %	0.01004 %	0.01229 %
0.10 °	0.00015 %	0.00011 %	0.00899 %	0.01350 %
0.30 °	0.00120 %	0.00105 %	0.00537 %	0.01887 %
0.95 °	0.01201 %	0.01062 %	0.00000 %	0.04193 %
3.00 °	0.12095 %	0.09733 %	0.05711 %	0.16350 %

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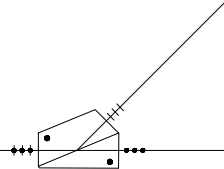The above table shows how reflectivity R_p of one surface is dependent upon small deviations δ from the Brewster angle α_B . The material Suprasil and a wavelength of 280 nm have been chosen for this example. The two columns to the right similarly show which reflections occur when this prism is used at 200 nm.

Bündelquerschnitt Beam cross-section	1 Stück Single	4 Stück Set of 4
10 x 10 mm	DBR 10	DBR 4.10
15 x 15 mm	DBR 15	DBR 4.15
20 x 20 mm	DBR 20	DBR 4.20

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Standard prisms are available in fused silica Suprasil 1; the ordering codes DBR 4.xx apply to the combination of 4 equal prisms often required in laser applications.

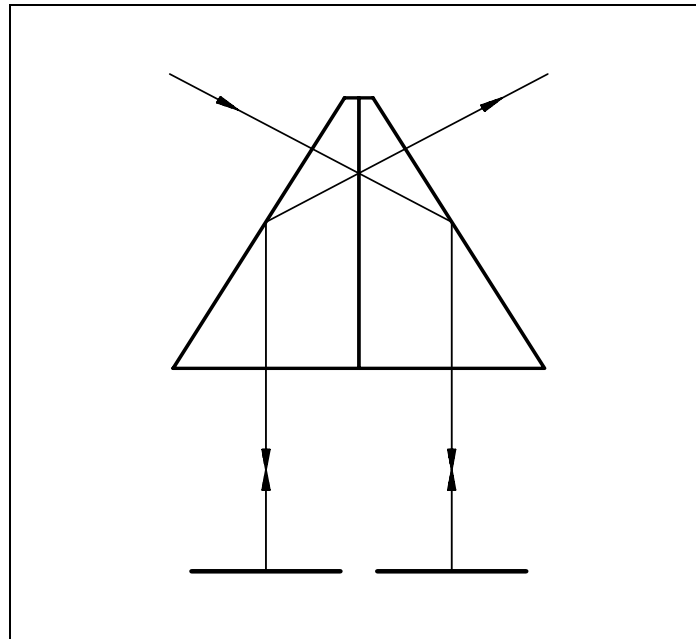
Ordering information:

When ordering, please enquire about prisms (wavelengths) which are currently available from stock.

Other sizes or materials on request

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Kösters Interference Double Prisms

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The prisms are made of fused silica Homosil and consist of two identical prisms halves which are cemented together. The flatness of the effective surfaces is $\lambda/10$. The angles of the prism halves are $30^\circ - 60^\circ - 90^\circ$, with angular accuracy $\pm 1'$ and angular equality $\pm 1''$. One long cathetus side is semi-transparent reflection coated with silver and a protective layer to $R = T = 45\%$ for the visible spectral range. The dimensions are listed in the table.

Dreiecksseitenlänge side length of triangle	große Kathete large cathetus	Fläche der kleinen Kathete surface of small cathetus	matte Höhe thickness	
20 mm	17 mm	10 x 10 mm	10	IKP 020
30 mm	26 mm	15 x 15 mm	15	IKP 030
40 mm	34 mm	20 x 20 mm	20	IKP 040
60 mm	52 mm	30 x 30 mm	30	IKP 060
100 mm	86 mm	50 x 50 mm	50	IKP 100

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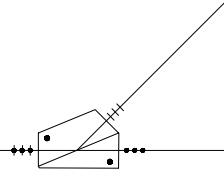
Available on special order:

Antireflection coatings of the entrance and exit faces on request

Beamsplitter coating with absorption-free dielectric layers on request


The two halves can also be supplied uncemented for use in special adjustable mounts. on request

Prism combinations for other interferometers on request

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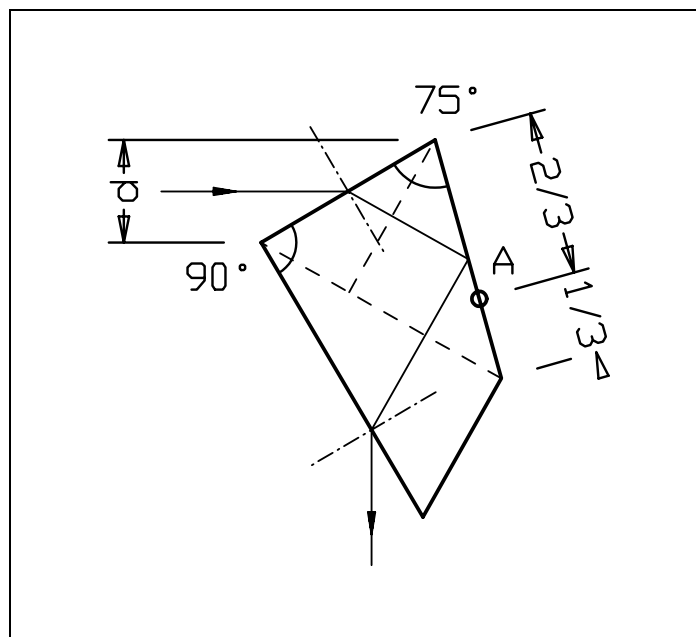These prisms are solid dispersing prisms with 90° deviation. They can be envisaged as a combination of two 30° dispersing prisms with one right angle reflection prism. The combination of the two 30° prisms works like a 60° dispersing prism and the 90° prism only does a pure reflection without dispersion.

The two 30° prisms (i.e. the 60° dispersing prism) are transmitted in the so-called „minimum of deviation“ by light of one wavelength. The entire prism deflects light of this wavelength by exactly 90° ; for this wavelength the spectroscopic angular dispersion is maximum.

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If the prism is rotated about axis A (see sketch), then a different wavelength is deviated by 90°. This way the prism can be tuned without the necessity to alter anything else in the geometry of the experimental setup.

The three polished surfaces of the prism produced from one block have a surface accuracy of $\lambda/10$. The angular errors are about 2'.

The following table lists a selection of Pellin-Broca prisms including individual beam cross-sections, material and reflection value R_s and R_p (per surface) of light polarized parallel or perpendicular to the plane of incidence (at some wavelengths λ).

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Other sizes and materials on request

Literatur:

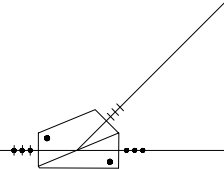
W. M. McClain, Appl. Opt. **12**, (1973), S. 153ff „How to Mount a Pellin-Broca Prism for Laser Work“

J.Reader, Appl. Opt. **12**, (1973), S. 1405, „Comment on: How to Mount a Pellin-Broca Prism for Laser Work“


Literature:

W. M. McClain, Appl. Opt. **12**, (1973), p. 153ff „How to Mount a Pellin-Broca Prism for Laser Work“

J.Reader, Appl. Opt. **12**, (1973), p. 1405, „Comment on: How to Mount a Pellin-Broca Prism for Laser Work“

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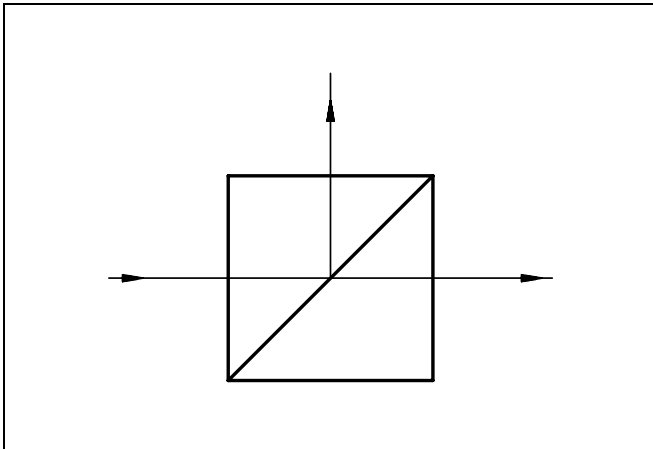The beamsplitter cubes are made of right angle prisms (crown glass BK 7) for a variety of spectral ranges.

The splitting layer is metallic (silver); the splitting ratio is 45:45% ($\pm 5\%$) for average polarization. The splitting layer slightly polarizes the splitted light: the transmission of the p-polarized component is about 5 to 15% higher than that of the s-component. The outer surfaces are broadband multilayer antireflection coated.

The surface accuracies of the right angle prisms are $\lambda/10$; the deformation of a transmitted wave is dependent of the homogeneity of the glass and often meets the quality of the surfaces. The prismatic deviation is not larger than about 1'.

Accessories:

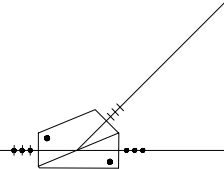
The cylindrical mounts holders describes on page 17 are suited for beamsplitters of sizes between 10 mm and 25 mm.

Available on special order:

Beamsplitter cubes with dielectric absorption-free splitting layer for one specified wavelength or a narrow wavelength range. on request


Beamsplitter cubes of fused silica for use in UV or in mid-range IR. on request

Beamsplitter cubes of raw glass with specified homogenities. on request

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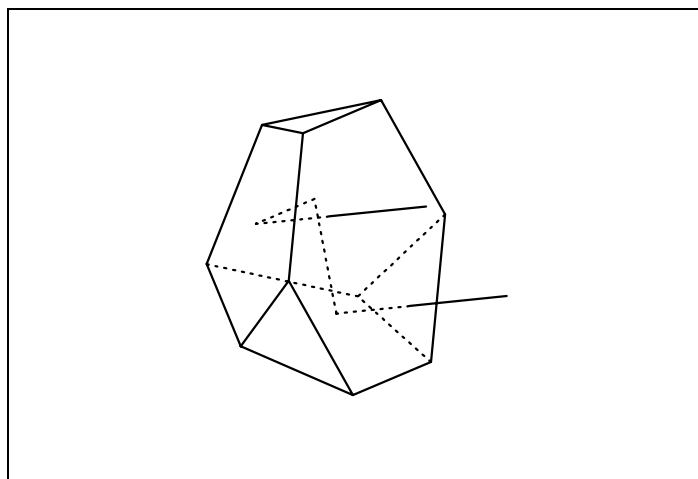The solid glass prisms are produced from fused silica Homosil or Suprasil and have sharp cube edges and surface flatness of $\lambda/10$. The deviation error is under 2". The shape of these prisms is either hexagonal or circular.

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Coatings:

Antireflection of the entrance faces with multilayer coatings for the visible. Minimal residual reflectivity about 0.2%STR 1

Antireflection of the entrance faces with double-layer coatings for a selected wavelength. Minimal residual reflectivity about 0.1%
STR 2

Reflecting surfaces coated with evaporated high-reflectivity metallic film to reduce the change of polarization
on request

Available on special order:

Prisms with other dimensions on request

Prisms with lower accuracies on request

The prisms can also be supplied mounted in a holder with or without a mounting post. on request

Hollow retroreflectors consisting of three flat mirrors of glass or Zerodur mounted by optical contact or cement. on request

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Known by the company name

Bernhard Halle Nachfl. GmbH

a team of scientists and high qualified fine optics engineers offers a large number of precision optical components. Especially the manufacturing of polarization optics is carried out with much experience. The present catalog contains our standard program. In addition to this program we offer special constructions which will be developed and realized according to your requirements together with our know-how.

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Optische Werkstätten *GmbH*