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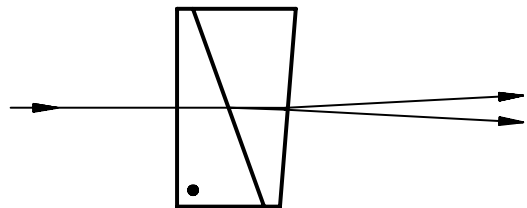
Hanle Wedge Depolarizers (Scrambler)

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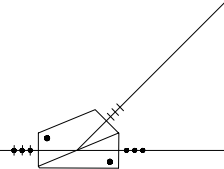
The depolarizer consists of two optically contacted wedges. At least one wedge is birefringent and affects the polarization state differently at different locations. The depolarization arises from the spatial superposition of rotated polarizations. The second wedge compensates for the prismatic deviation (as good as possible). The divergence of the two emerging beams polarized at right angles to each other is about 4 arc minutes. Flatness of the outer surfaces is $\lambda/10$. The total thickness is about 6 to 10 mm.

The depolarizer influences the polarization of light sources or of a device's own polarization so that the average degree of polarization is close to zero. Thus for each respective polarization the depolarizer must be adjusted to the most favourable orientation achievable by experimentation. (Usually it is a good starting value to orient the divergence under 45° to the prevailing polarization direction of the incident light) The effectiveness of the depolarizer increases with the size of the beam cross-section.

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Type I:

Crystal wedge of synthetic quartz, compensator wedge of fused silica Suprasil. Spectral range about 180 nm - 2.5 μ m. unmounted.

Querschnitt (achteckig) cross-section (octogonal)	synthetischer Quarz synthetic quartz
10 x 10 mm	PDH 10
15 x 15 mm	PDH 15
20 x 20 mm	PDH 20
25 x 25 mm	PDH 25

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Type II:

Crystal and compensator wedge of MgF₂, crystal axes crossed, spectral range about 125 nm - 7 μ m, unmounted

Cross section 15x15 PDH 15.00

The depolarizers can be supplied in a mount with a mounting post. Mounted in such a holder the depolarizer can be rotated 360° and locked into any position. The clear diameters of the holders are 9 mm, 14 mm, 19 mm and 24 mm; the outer diameters are 30 mm, 30 mm, 36 mm and 38 mm. PDH 0

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Literature:

translated title: „Measurement of the degree of polarization of spectral lines.“